



Mast Group Ltd. Mast House, Derby Road, Bootle, Merseyside, L20 1EA United Kingdom

Tel: + 44 (0) 151 472 1444 Fax: + 44 (0) 151 944 1332 email: sales@mast-group.com Web: www.mast-group.com



Mast Diagnostica GmbH
Feldstrasse 20
DE-23858 Reinfeld

Tel: + 49 (0) 4533 2007 0 Fax: + 49 (0) 4533 2007 68 email: mast@mast-diagnostica.de Web: www.mast-group.com

Germany

Mast Diagnostic

12 rue Jean-Jacques Mention CS91106, 80011 Amiens, CEDEX 1 France

Tél: + 33 (0) 3 22 80 80 67 Fax: + 33 (0) 3 22 80 99 22 email: info@mast-diagnostic.fr Web: www.mast-group.com



# **Blood Agar Base**

### **DM100**

## **Intended Use**

A general purpose medium that gives typical haemolytic reactions when the medium is supplemented with sterile defibrinated blood.

### **Contents**

See pack label.

### Formulation\*

Material:	Concentration in medium:
Peptone mixture	16.0g/litre
Yeast extract	2.0g/litre
D-glucose	0.5g/litre
NaCl	7.0g/litre
Agar	12.0g/litre
Final pH: $7.3 \pm 0.2$	

## Storage and shelf life

All dehydrated culture media containers should be kept tightly closed and stored in a dry place at 10 to 25°C until the expiry date shown on the pack label.

#### **Precautions**

For *in vitro* diagnostic use only. Observe approved hazard precautions and aseptic techniques. To be used only by adequately trained and qualified laboratory personnel. Sterilise all biohazard waste before disposal. Refer to Product Safety Data sheet (available on request or via MAST® website).

### Materials required but not provided

Standard microbiological supplies and equipment such as loops, MAST® selective supplements, swabs, applicator sticks, incinerators and incubators, etc., as well as serological and biochemical reagents and additives such as blood.

## **Procedure**

- Refer to pack label for quantities and volumes required. Prepare MAST® Blood Agar Base (DM100D) by suspending the powder in distilled or deionised water. For sachet packs, dissolve the entire contents of the sachet in the volume shown on the label.
- 2. Autoclave at 121°C (15 p.s.i.) for 15 minutes.
- Cool to 50 to 55°C and hold at this temperature in a water bath. Add 5 to 7% sterile defibrinated horse or sheep blood where required. Heated blood agar (chocolate) can also be prepared. Mix well before pouring.
- 4. Pour culture plates (15 to 20ml per plate) and allow to

- 5. Prepared culture plates may be used immediately or stored in plastic bags at 2 to 8°C for up to one week before use.
- 6. Inoculate plates by surface plating, streaking out for single colonies.
- Incubate plates aerobically for 18 to 24 hours and anaerobically for up to 72 hours at 35 to 37°C (or alternative temperatures according to the methodology followed).

## Interpretation of results

After incubation record growth of organisms. Typical characteristics to note include colony size and morphology, pigmentation and haemolysis.

### Quality control

Check for signs of deterioration. Quality control must be performed with at least one organism to demonstrate expected performance. Do not use the product if the result with the control organism is incorrect. The list below illustrates a range of performance control strains which the end user can easily obtain.

Test Organisms	Result
Escherichia coli	Growth
ATCC® 25922	
Staphylococcus aureus	Growth,
ATCC® 25923	β-haemolysis
Streptococcus pneumoniae	Growth,
ATCC® 6305	$\alpha$ -haemolysis
Streptococcus pyogenes	Growth,
ATCC® 19615	β-haemolysis

### References

Bibliography available on request.